MILK DOCTOR BETA

Kit Compositions

Test Strips: 100 pieces Sample cup: 100 pieces Dropper: 100 pieces

TEST PROCEDURES



1

Take the milk sample with a Dropper up to the level of the indicated line, and dispense all amount of the milk into the plastic Cup.

Do this operation Twice.





2

Jiggle the Plastic Cup lightly to spread the sample on the bottom surface of plastic cup







3

Soak the Dipping part of Test Strip in the milk sample of the plastic Cup.

* Do not place the strip vertically







MILK DOCTOR BETA

TEST RESULT



Think Quality & Create Value